

M-MLV RTase (M-MLV Reverse Transcriptase)

Cat no. LDG0006RF

Product Overview

Package component

Specification	Item	Amount
20,000 U	M-MLV Reverse Transcriptase	1 vial (200 U/μL)
	5X M-MLV Reverse Transcriptase Reaction Buffer	1 vial (1 mL)
50,000 U	M-MLV Reverse Transcriptase	1 vial (200 U/μL)
	5X M-MLV Reverse Transcriptase Reaction Buffer	1 vial (1 mL)

Description

Moloney Murine Leukemia Virus (M-MLV) Reverse Transcriptase is an RNA-dependent DNA polymerase that synthesizes the first strand of complementary cDNA from a single-stranded RNA template with hybridized primer. This kit features high activity formulation of M-MLV RT and 5X reverse transcription buffer which are capable of full-length cDNA synthesis and high cDNA yields.

Source

Escherichia coli

Activity

One unit of M-MLV Reverse Transcriptase is defined as the amount of the enzyme incorporates 1 nmol of dTTP into acid-insoluble product in 10 minutes at 37°C.

Storage buffer

The enzyme is supplied in 20 mM Tris-HCl (pH 7.5), 200 mM NaCl, 0.25 mM EDTA, 0.01% NP-40 (v/v), 2.5 mM DTT, 50% glycerol (v/v).

5X M-MLV Reverse Transcriptase Buffer

250 mM Tris-HCl (pH 8.3), 15 mM MgCl₂, 375 mM KCl, and 50 mM DTT.

Storage and Stability

Stored at -20°C. Avoid repeated freeze/thaw cycles.

Procedure

First strand cDNA synthesis:

- Place all required reagents to a nuclease-free microcentrifuge tube and following the order suggested below.

Component	Amount	Final concentration
Oligo (dT) 12-18 (50 μM) or random primer mix (60 μM)	1 μL	-
Total RNA template	X μL	1 μg
Nuclease-Free H ₂ O	Y μL	-
Total reaction volume	10 μL	-

- Heat the tube to 65°C for 10 minutes to denature the secondary structure within RNA template. Immediately cool the tube on ice for 1 minute and centrifuge briefly in microcentrifuge. Add the following components to the annealed primer/RNA template, prepare on ice.

Component	Amount	Final concentration
5X Reverse Transcriptase Reaction Buffer	4 μL	1X
10 mM dNTPs mix	1 μL	0.5 mM each
RNase Inhibitor	X μL	20 U/rxn
M-MLV RTase	1 μL	200 U/rxn
Nuclease-Free H ₂ O	Y μL	-
Total reaction volume	20 L	

- Incubate at 37°C for 1 hour. The extension temperature may be adjusted from 37°C to 42°C.

4. Inactivate the reaction at 65°C for 20 minutes. The cDNA products should be store at -20°C.
5. Reaction preparations may be scaled up or down proportionately.

Important notes

- (1) After the reaction is complete, M-MLV RTase can be inactivated by incubation at 65°C for 20 minutes.

Disclaimer

This product is for research use only and is not intended for diagnostic use.

LEADGENE BIOMEDICAL, INC.

No.9, Ln. 147, Zhengbei 1st Rd., Yongkang Dist., Tainan City 710, Taiwan R.O.C. TEL: +886-6-2536677 FAX: +886-6-2531536
www.leadgenebio.com