



Product Information & Manual

Information of other products is available at: www.leadgenebio.com

Leadgene® StablePlus™ 2X Colorimetric RT-LAMP Master Mix

Cat no. LDG0024RF

Product Overview

Package component

Item	Content
StablePlus™ 2X Colorimetric RT-LAMP Master Mix	1 vial (1 mL)

Description

Leadgene® StablePlus™ 2X Colorimetric RT-LAMP Master Mix is an optimized master mix for reverse-transcription loop-mediated isothermal amplification (RT-LAMP) reactions. This product is a dual enzyme system, providing a rapid and sensitive detection in one pot.

The amplified products can be visually observed by a metal indicator for calcium, hydroxynaphthol blue (HNB). Positive reactions will appear a color change from violet to sky blue, while negative reactions won't.

The StablePlus™ version contains nucleic acid stabilizing agent to protect the amplified products.

Storage and Stability

Stored at -20°C. Avoid repeated freeze/thaw cycles.

Procedure

The following procedure is a general guideline for RT-LAMP reaction. To maintain an RNase-free environment, always wear disposable gloves, and use laboratory consumables and water of nuclease-free grade during the whole experiment course.

RT-LAMP reaction set-up:

- 10X LAMP primer mix

Component	10X concentration	Final concentration
FIP	16 µM	1.6 µM

BIP	16 µM	1.6 µM
F3	2 µM	0.2 µM
B3	2 µM	0.2 µM
LOOP F	8 µM	0.8 µM
LOOP B	8 µM	0.8 µM

- An overview of the reaction set-up is listed in the table below. Place all required reagents **on ice**. Distribute appropriate volumes into each tube before adding template.

Component	Amount	Final concentration
StablePlus™ 2X Colorimetric RT-LAMP Master Mix	12.5 µL	1X
10X LAMP primer mix	2.5 µL	1X
RNA template	1-2 µL	variable
Nuclease-Free H ₂ O	X µL	-
Total reaction volume	25 µL	-

- Add target RNA template to the detection tube. Gently mix the reaction thoroughly to achieve uniform distribution and avoid making bubbles.
- Incubate at 65°C for 30-60 min.
- After LAMP reaction complete, the enzyme can be inactivated by heating at 80°C for 10 min.

Important notes

Primer concentration

Primer concentration can be titrated between 0.25X – 1X if undesired background signal appeared.

Disclaimer

This product is for research use only and is not intended for diagnostic use.

LEADGENE BIOMEDICAL, INC.

No.9, Ln. 147, Zhengbei 1st Rd., Yongkang Dist., Tainan City 710, Taiwan R.O.C. TEL: +886-6-2536677 FAX: +886-6-2531536
www.leadgenebio.com